


Function Button Press

ON/OFF	Power up		1 x long
	Power down		3 x short
	Activate key lock		1 x long
	Deactivate key lock		

Menu operation




The active mode is highlighted by a ✓

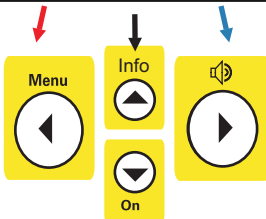

Selection of a digit, decimal separator or prefix: ← →


Changing its value: ↑ ↓

Ratemeter Mode *(Fast scanning for surface contamination)*

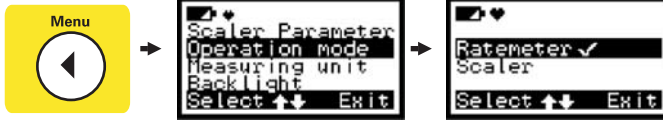

1 Battery Level

3 Bar graph display of the count rate in logarithmic scale

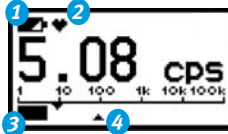

2 Icon indicates, that the unit is processing data

4 Alarm threshold setting:
Level 1: ▼ Level 2: ▲

5 Quick count rate change.
Wait until the count rate is stable
Increasing count rate ↗
Decreasing count rate ↘



6 Count rate indication by sound (pulse)

2x




Scaler Mode *(Highly sensitive measurements of prepared samples)*

The RadEye B20 / B20-ER can be set to display the net count rate by automatically subtracting the stored background data. See the manual for details.



1 Preset measurement time


2 Press  to start the measurement

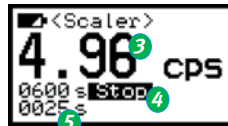

3 Average value based on the elapsed time

4 Early termination of the measurement

5 Elapsed measurement time

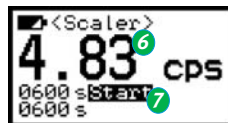

6 Measurement result

7 Press  to start the next measurement